

PANalytical X-Ray Diffractometer Standard Operating Procedure

Version 2.1 (Corrected on March 2024)



Training Rules:

First time: watch trainer and take notes

Second time: do by yourself with supervision following your notes

Third time: do by yourself with supervision without help (Dr. Bykov must approve the authorization for independent use)

After the third time, you are completely independent and responsible for the instrument.

Operating Rules:

1. Only authorized people can use XRD.
2. You must sign up for time on the XRD using the online system:
<http://corelabs.ccnycuny.edu/>
3. All samples must be recorded in XRD logbook including the date, time, user, sample name, program name, and filename.
4. If you have any problems, please DO NOT troubleshoot yourself - Call:
Alexey Bykov 212-650-5548 (Laboratory) or 646-725-0270 (cell phone)

Outline:

1. Turn on the instrument
 2. Tube breeding
 3. Make a new program
 4. Typical operation with fixed holder
 5. Typical operation with spinner holder
 6. Export data
 7. Turn off the instrument
-

1. Turn on the Instrument:


1.1 Press power switch on the water chiller.

1.2 If PC is ON, double click on **X'Pert Data Collector** icon.




If software is already running but generator status is OFF (Tension = 0 kV and Current = 0 mA) skip steps 1.2-1.5 and go to step 1.6. If generator status is also ON (Tension > 20 kV, Current > 5 mA) go to step 2.3.


1.3 Type user name: **User-1** and password: **galaxy** in Login dialogue box.

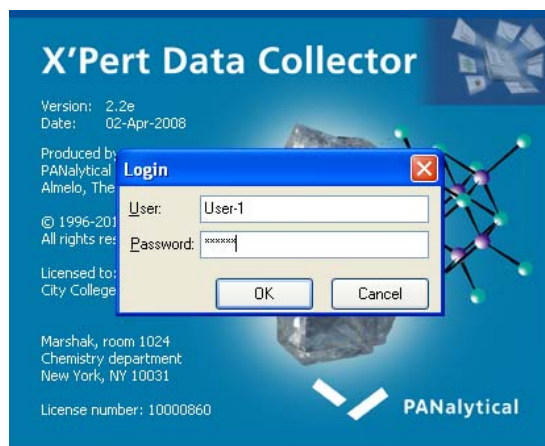


1.4 Go to **Instrument** menu and click **Connect**.

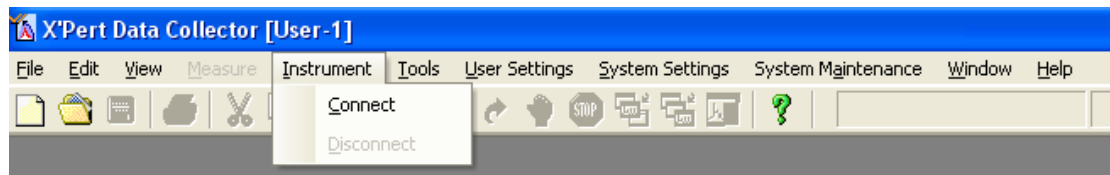

1.5 Select **Stage Flat Sample** for regular measurements or **Sample Spinner** if you need to use sample changer and/or rotate your sample during scanning procedure and press **OK** (To switch sample holders go to **Tools** menu → **Exchange Sample Stage** with following connection procedure

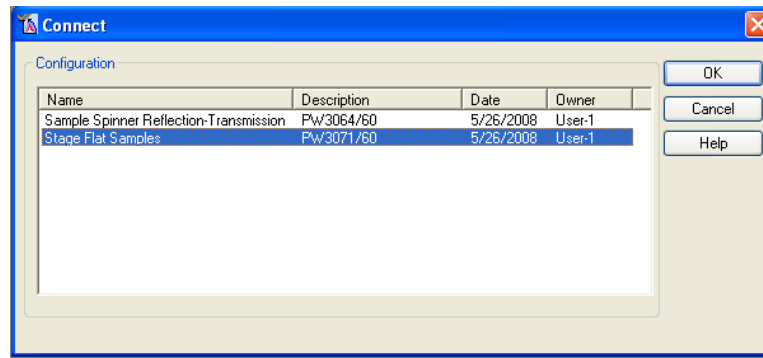


1.6 Double click on **X-ray**. Check the **Generator on** box at the **Instrument Setting**, press **Apply**, and wait until tube power reaches 30 kV and 10 mA .

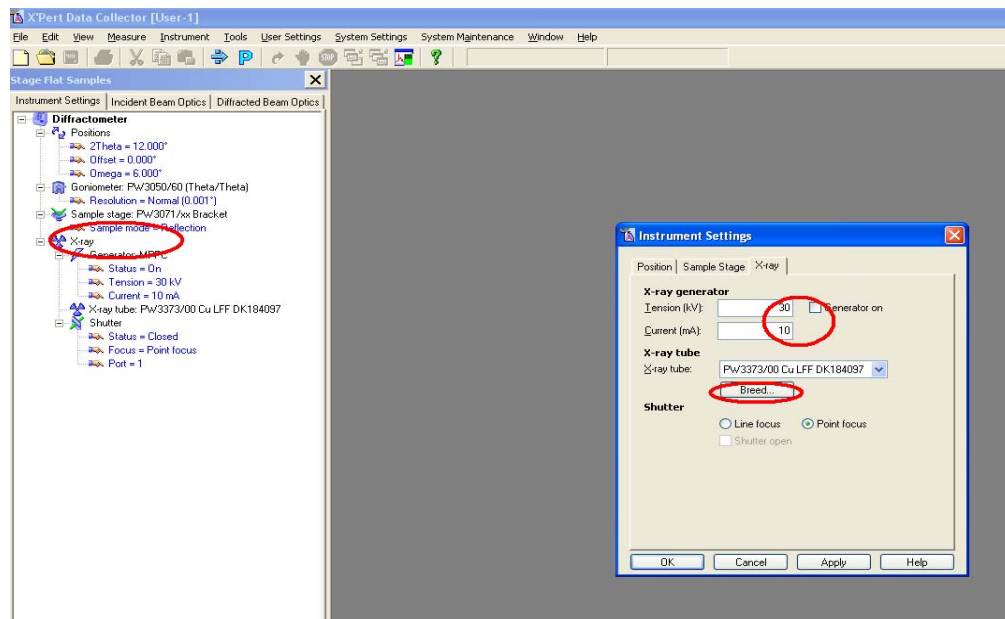

2. Tube Breeding

If X-rays have been turned off for an extended period of time, tube breeding (conditioning) is necessary. Check the log book first to see when last time the instrument was used. Breed at normal speed if the instrument was left idle for 100 hours or more (breed takes about 30 minutes). Breed at fast speed if the instrument was used recently (within the last 100 hours). In this case, tube breeding takes about 6 min.

2.1 Click **Breed**. Choose **fast** or **normal** in **Tube Breeding** box and mark **OK**

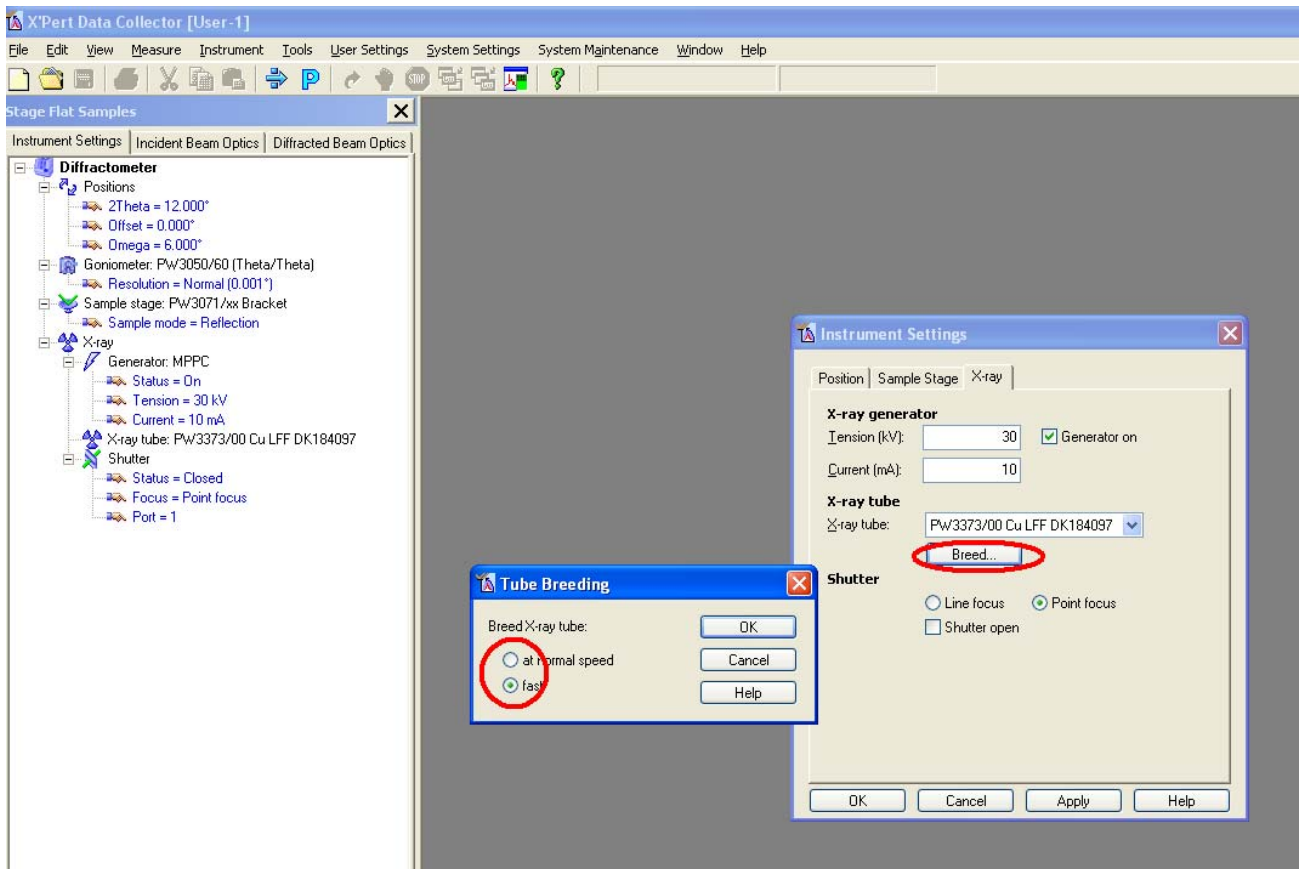


2.2 Tension during breeding procedure goes up to 60 kV then down to 40 kV.

After breeding the parameters of the generator will be **40 kV and 10 mA**

2.3 On the **Instrument Settings** window set measurement parameters **40 kV** and **40 mA**. Click **Apply** then **OK**.

3. To Make a New Program:

3.1 Go to **File** menu and click **Newprogram**.

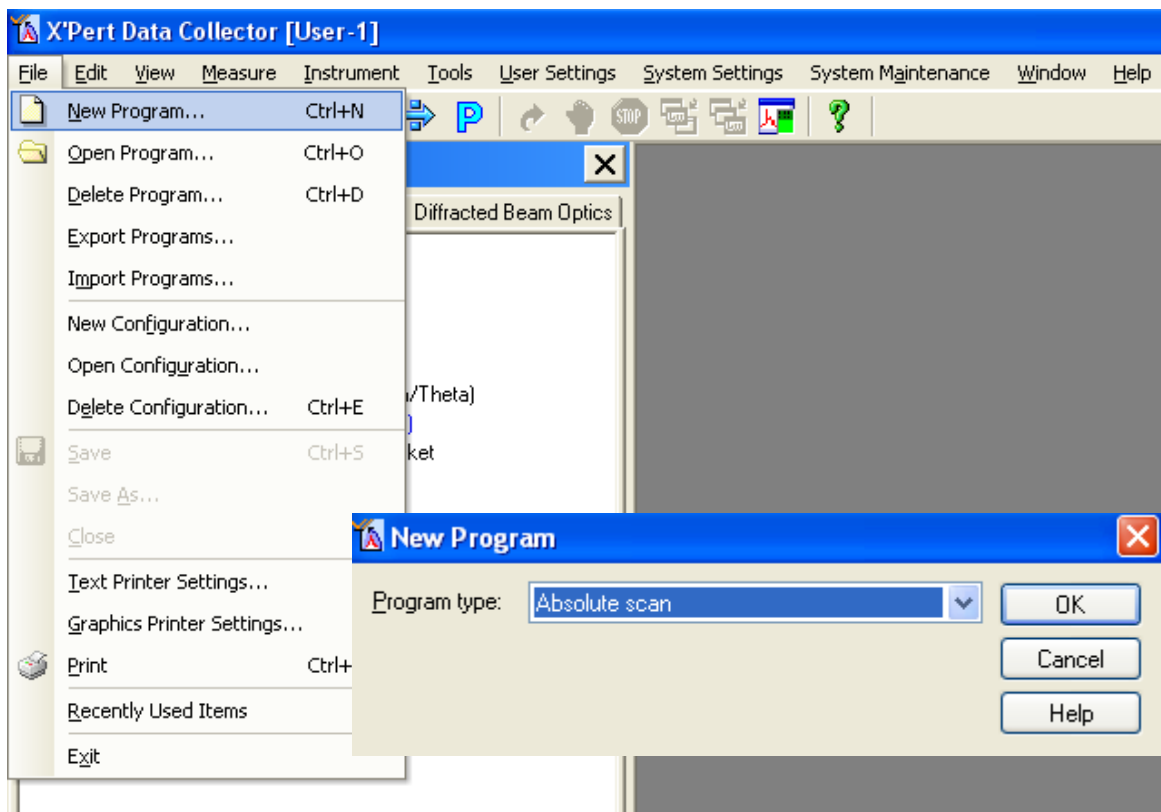


Click **OK** for **Absolute scan** program type.



If you have already created your program in previous sessions, use **Open Program** in **File** menu and make parameters changes if you need it, then go to step 4 (skip step 3.2 -

3.3)

3.2 Choose your scanning parameters on the Prepare absolute scan window.

- In Configuration, select stageflatsamples (for fixed holder) or Sample reflection Transmission (for spinner holder)
- In Scan Axis, select the type you want, a typical scan for powders is Gonio.
- Click on Repetition, select then Single scan or other scan type.
- Click Settings
- Click Detector: Actual, choose Pixel(2)(window shows the parameters of PIXcel detector: *Active length* (angle) and *Number of active channels*)
- Click Apply, then OK

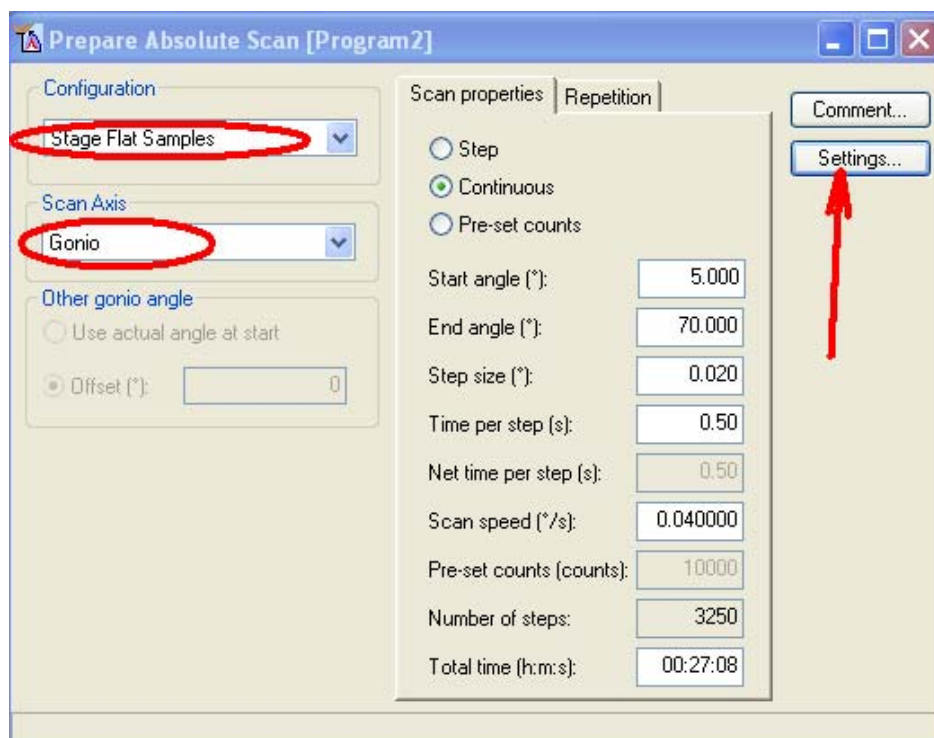

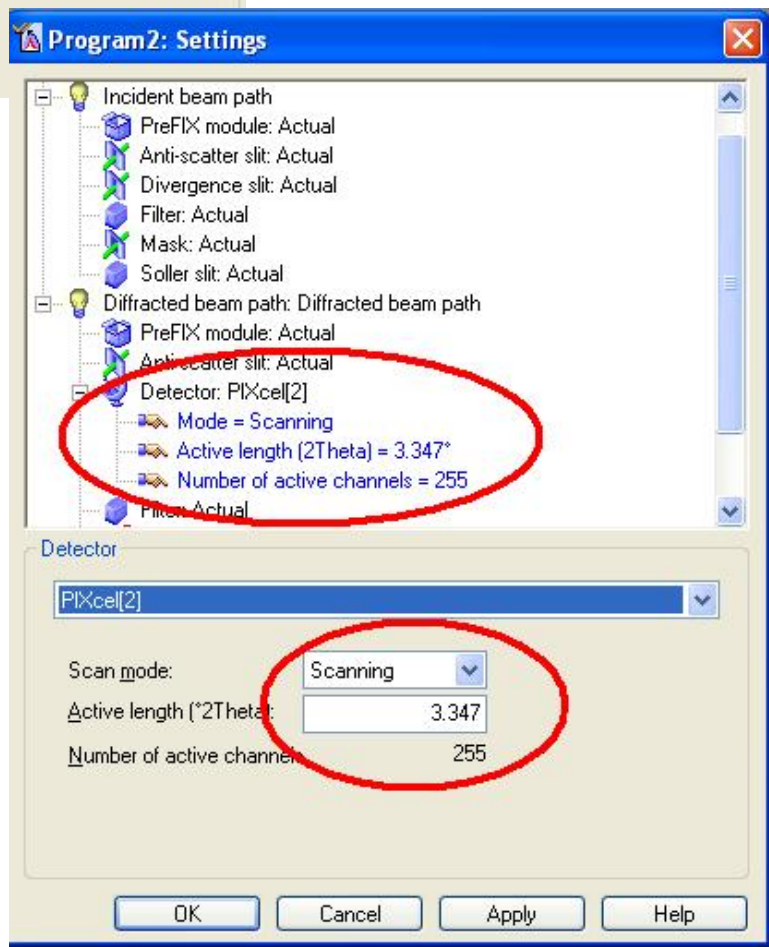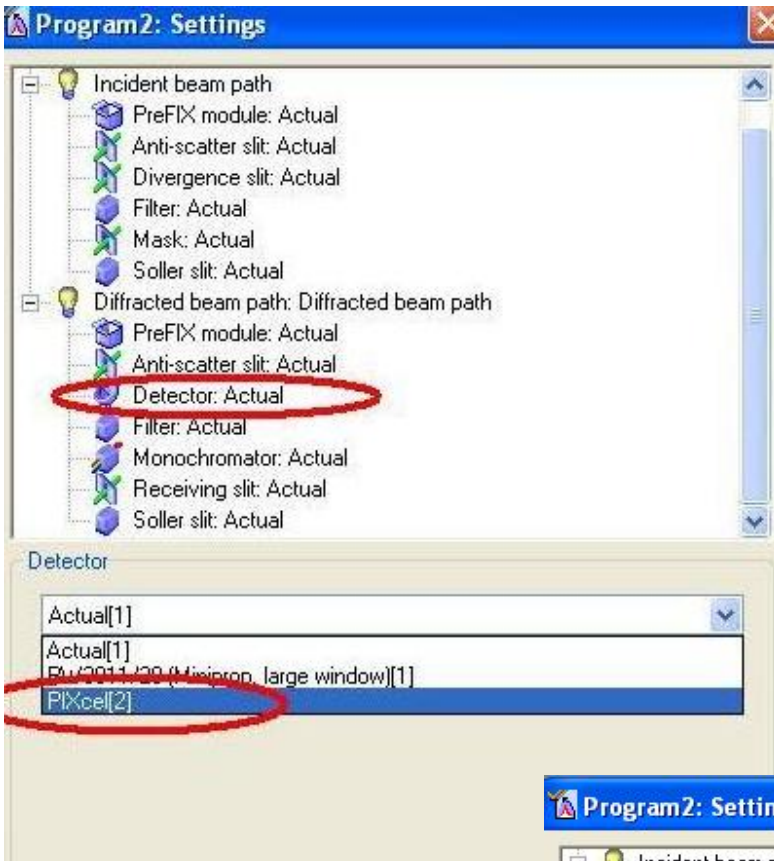

PIXcel detector



PIXcel detector has been developed by PANalytical as an advanced two-dimensional photon counting device incorporating the latest in solid-state pixel technology. In our system, PIXcel detector works in 1D (one dimensional) mode. The detector simultaneously counts photons in 2Theta range = 3.347° by 255 active channels. This is a maximum detector's active length. The active length may be decreased down to 0.013° and only one channel will be active. At this setting the detector is converted to 0D mode and works like point detector.

- Click Scan properties tab to edit the scanning parameters. Only Continuous scan is allowed by PIXcel detector. Set your angle range, step size, time per step (or scan speed). Based on our tests of standard samples we recommend for regular measurements step size ~ 0.05 - 0.01° and scan speed ~ 0.1 - $0.2^\circ/\text{sec}$. The total scan time usually should not exceed 10 min. (5 min is optimal total time for 10° - 60° Two Theta range)

PIXcel detector at a maximum value of active length (2Theta range = 3.347°) allows step size range from $\sim 0.002^\circ$ to 0.512° and scan speed up to $1.27^\circ/\text{sec}$

3.3 Save experiment

- Go to main menu, select File, click Save as or click close the Prepare Absolute Scan window and press Yes to save the file (use your first name or any other title (Ex: Mineral_2) and then click OK).

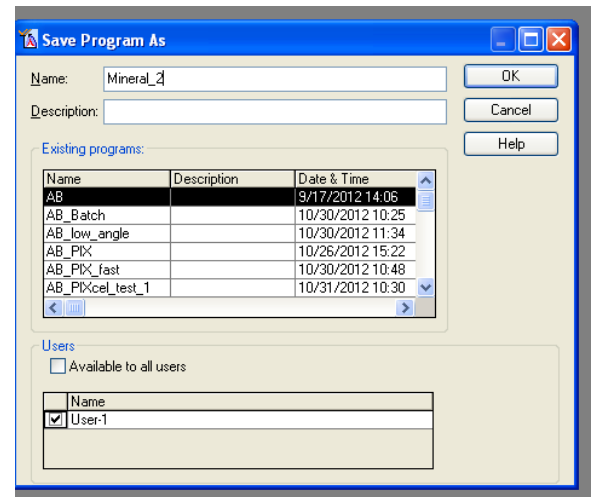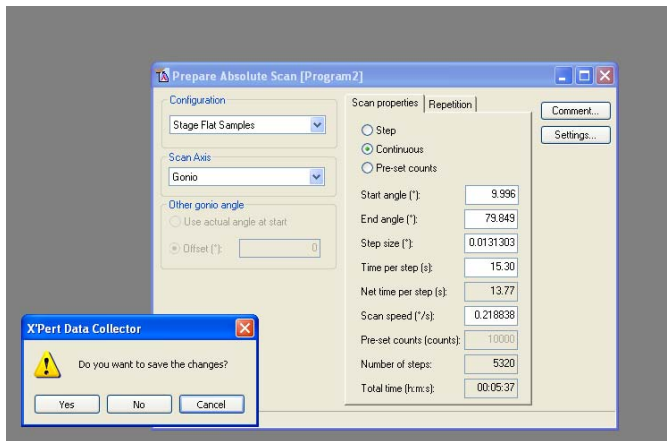



For previous saved program with running XRD you may **start here**.

4. Typical Operation with Fixed Holder

4.1. Load XRD sample slides on the fixed sample holder.

4.2. Go to main menu, click **File**, choose **Open program** and check the program settings.

- 4.3. Go to main menu, Select **Measure**, click **Program**, and **OK** after you choose your program.
- 4.4. Click the folder icon on **Start** window (If you are a new user please create a folder under your group folder in **XPert Data** directory)
- 4.5. Type in the file name, then click **Save**
- 4.6. Click **OK** on **Start** window. The XRD instrument starts scanning.

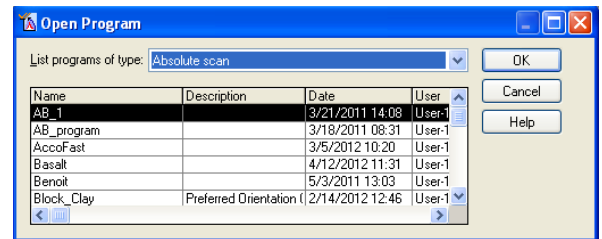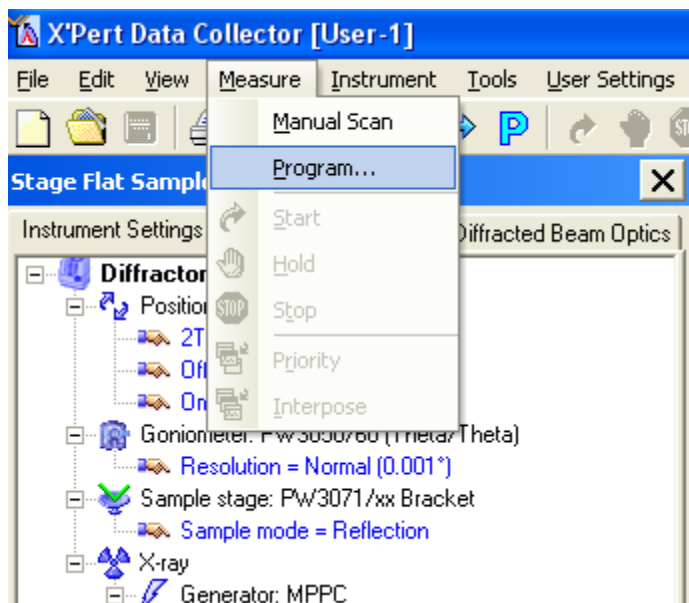

Start

Program Name: AB_1
 Type: Absolute scan
 Description:

File Name:
 Folder: C:\XPert Data
 Comment:

Sample ID:
 Name:
 Prepared by:

Position
 Diffractometer
 2Theta (°): 5.000 Phi (°): X (mm):
 Offset (°): 0.000 Psi (°): Y (mm):
 Omega (°): 2.500 Z (mm):
 Reflection Unit cell: h k l:

OK Cancel Help

Start


Program Name: AB_1
 Type: Absolute scan

Select Filename

Save in: X:\Pert Data






File name: cement
 Save as type: Measurement Files (*.xrdml)

OK Cancel Help

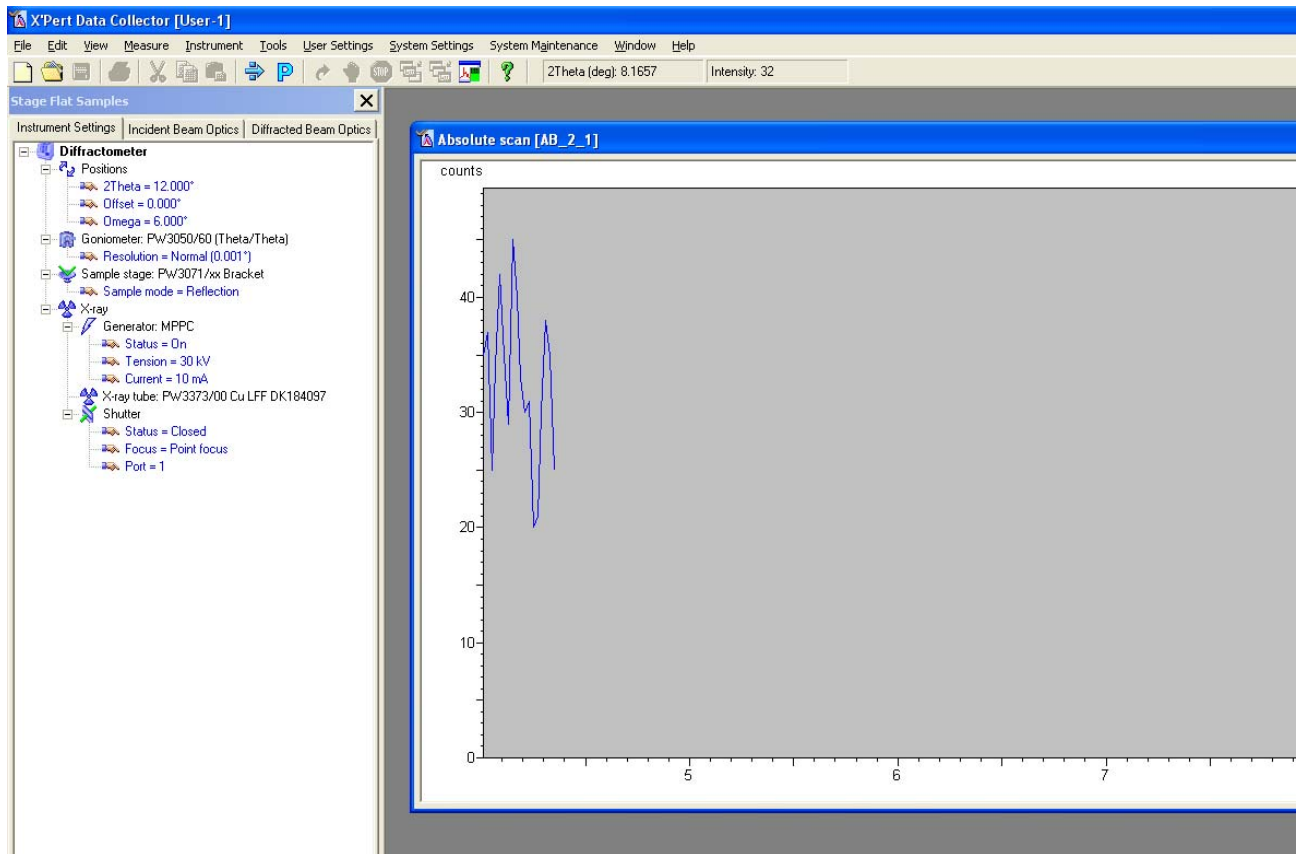


5. Typical Operation with Spinner Holder



5.1. To switch sample holders you should disconnect XRD **Instrument – Disconnect**, then go to **Tools** menu → **Exchange Sample Stage**, choose **Reflection Transmission Spinner** at **Change Sample Stage** window, press **Apply**. Follow the instructions of the **X'Pert Wizard**

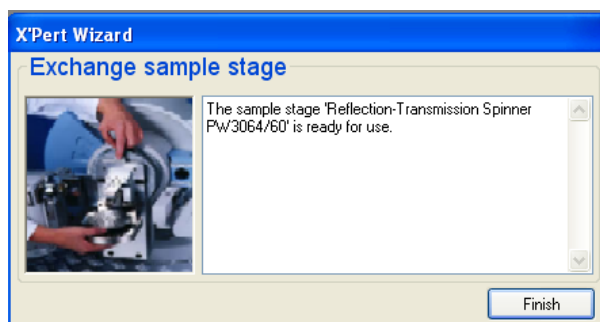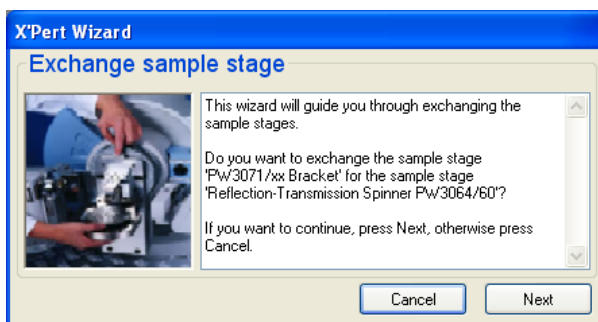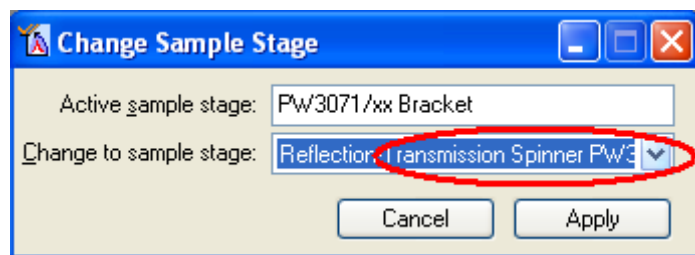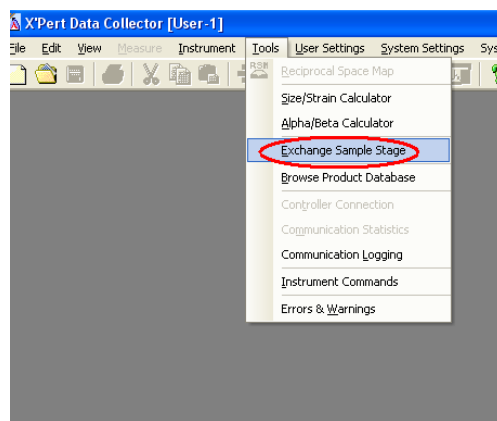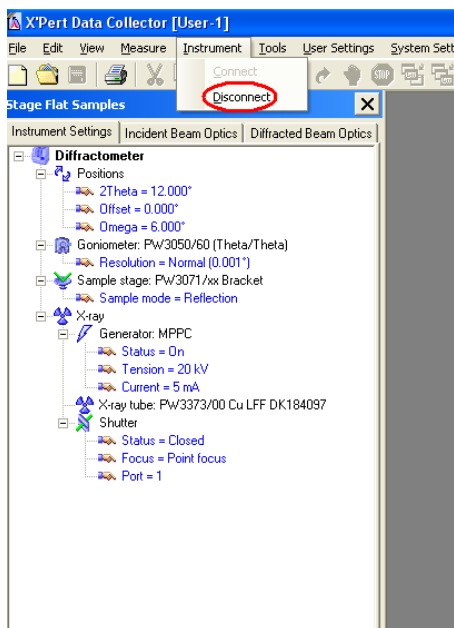


The Fixed Stage is mounted to the base by four screws. Unscrew all of them and carefully remove the stage. Take the Spinner Holder and attach to the base. Mount the holder using three screws only.

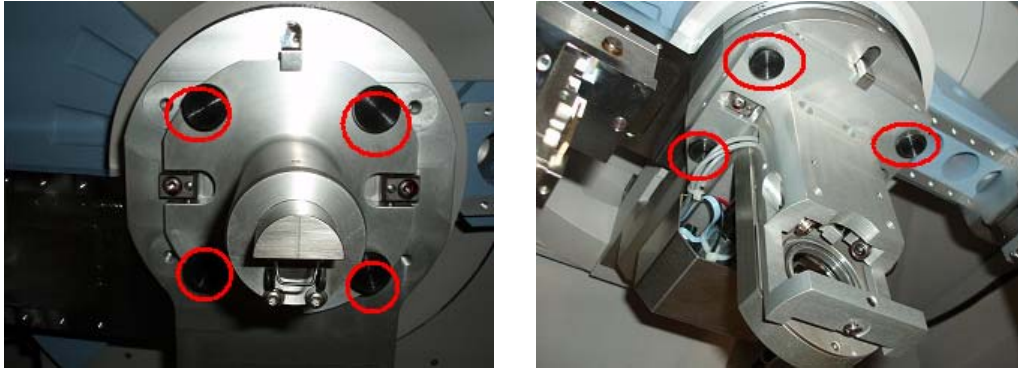

- Connect the instrument (Go to Instrument menu and click Connect).
- Select Sample stage Reflection Transmission in the Connect window and click OK

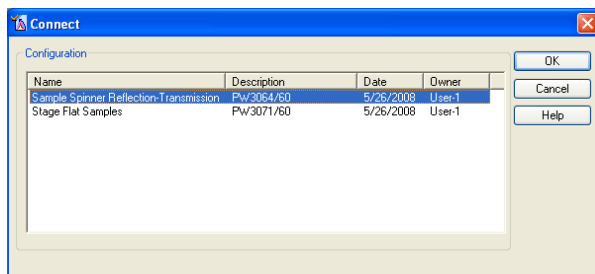
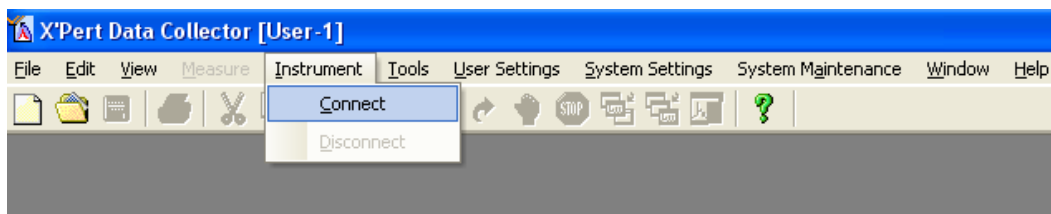


If the Sample Changer is on corner position it's status will be Unavailable

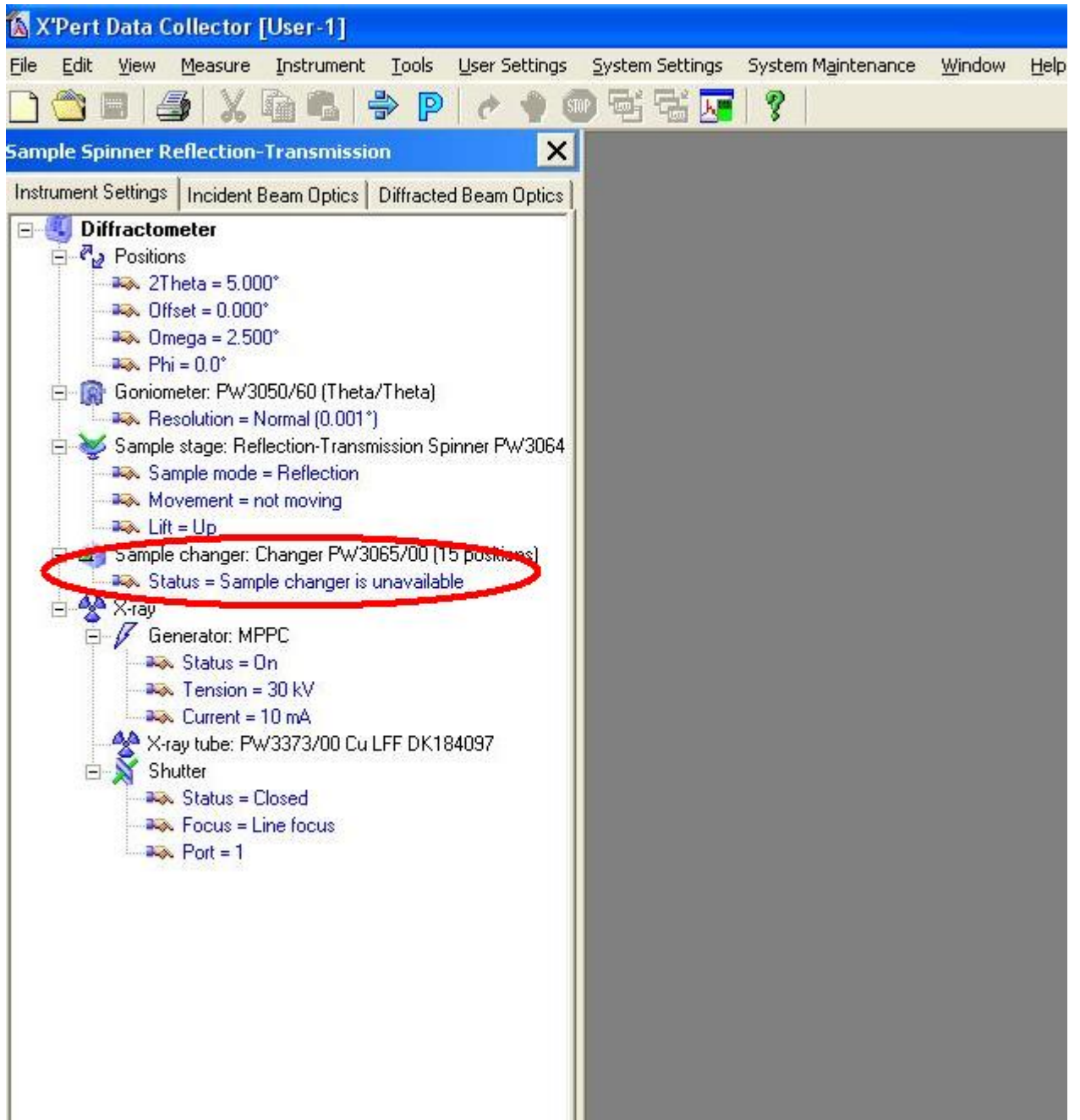


- Move Sample Changer from corner to center position

This will be done by XRD facility manager only !!!

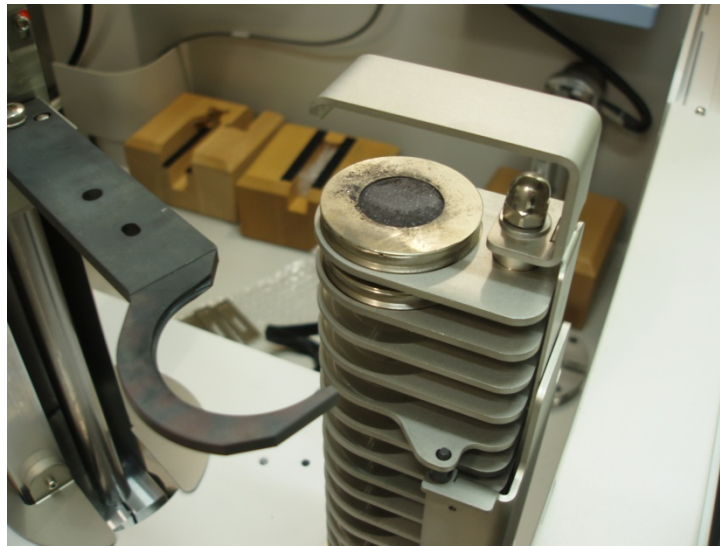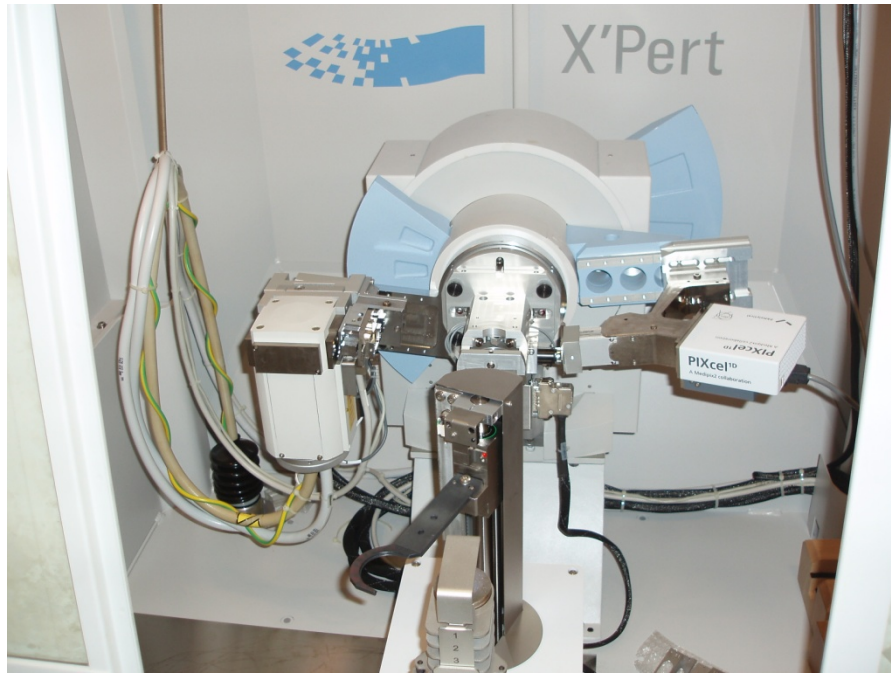

When Sample Changer will be fixed in center position it's status will be changed from Unavailable to No sample loaded. If there is no changes in Sample Changer status occurred, click on Sample Changer to open Instrument setting window, mark Sample changer available box and press Apply.

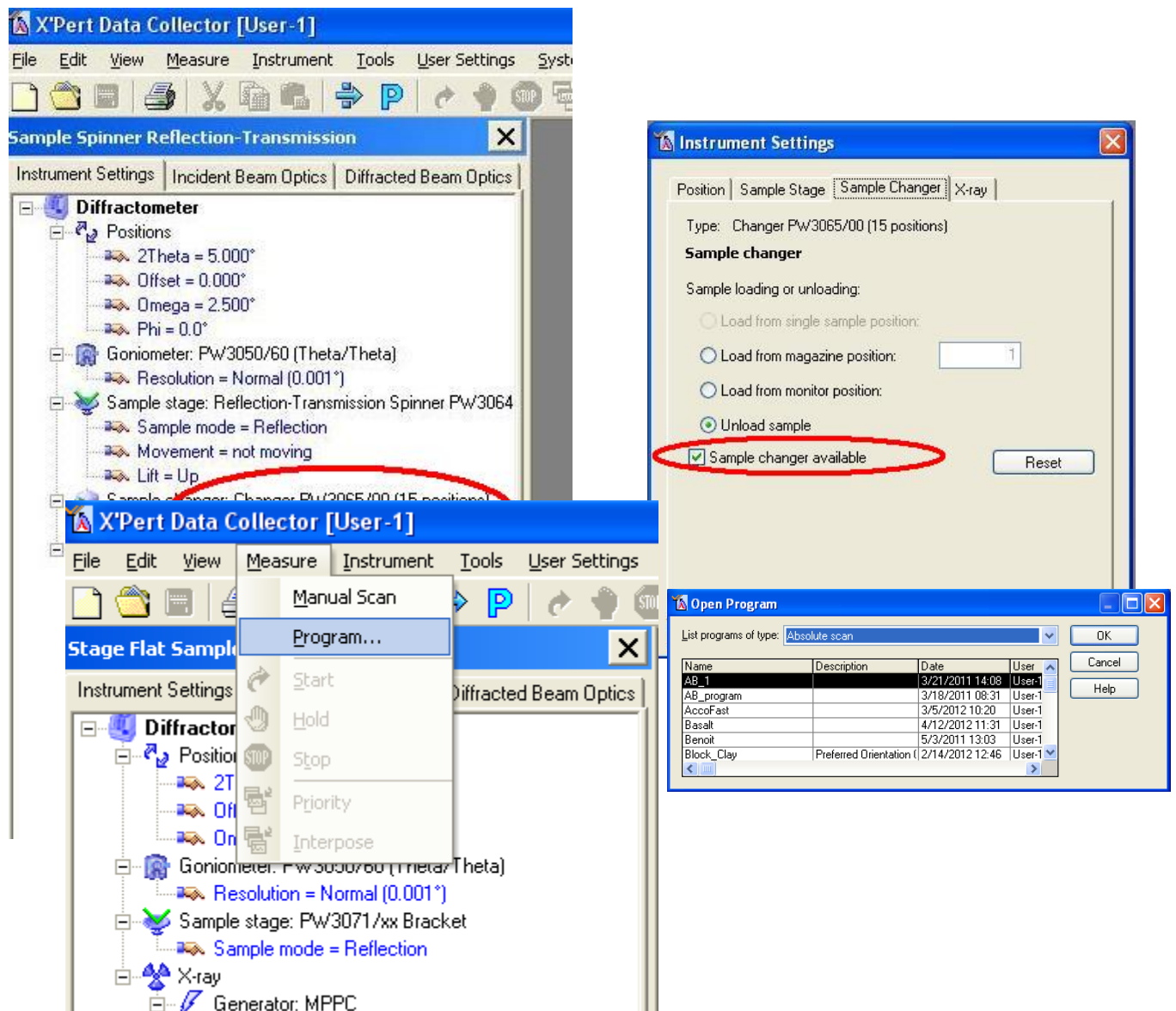

5.1. Load your sample on position 1 (top) of the Sample Changer.

5.2. Go to the **File** menu then choose **Open program**. The program settings are appropriate to **samplespinnerReflection-Transmission** mode (**Configuration** – **Sample Spinner Reflection**, **Scan Axis** – **Gonid**. Click **OK**

5.3. Select **Measure** and click on **Program**, then **OK** after you choose your program.

5.4. Click the folder icon on **Start** window (If you are a new user please create a folder under your group folder in **XPert Data** directory)

5.5. Type in the file name, then click **Save**

5.6. Click **OK** on **Start** window. The sample will be loaded in spinner and XRD instrument starts scanning. After completing measurements the Changer will unload spinner holder placing your sample back to position

6. To export data:

6.1 On the desktop, double click on **High Score Plus** icon to open analyzing software.

6.2 Go to **File** menu then **Open** select the file from your folder.

6.3 In the **ListPanel**, select **Anchor Scan Data** tab, mouse right click while marker on the excel chart. To get XY data, select **Save list as**, and name your file. Save the data in .CSV or .txt format file.

6.4 To export data in different format, go to **File** menu and select **Save As** and select the desired data format, such as: RD, UDF, DAT...etc.

Data saved in C:\X'Pert Data\ automatically

7. To Turn off the Instrument:

7.1 Double click on **X-ray** on **Stage Flat Sample** or **Sample SpinnersStageReflection-Transmission** on left side of **Data Collector** window

7.2 Change tube power to 20 kV and 5 mA in **Instrument Settings** box.

7.3 Click then wait for ~1-2 min, uncheck the *Generator* box and click .

(Do not turn off the key on instrument panel).

7.4 Turn off the cooling water switch on chiller (recirculator).